# DNBSEQ™SERVICE OVERVIEW **Metagenomic Sequencing**



# **Service Description**

Metagenomics studies of genetic material that is directly recovered from environmental samples, have benefitted greatly from advanced NGS technology as a method for the exploration of microbial biodiversity. BGI's metagenomics survey service applies whole genome shotgun sequencing of DNA isolated from environmental samples, with the advantages of high throughput and high coverage. It can provide information not only on species composition and abundance, but also on functional genes, gene differences between samples, metabolic pathways and gene resource mining for bioactive products.

# **Project Workflow**

We care for your samples from the start through to the result reporting. Highly experienced laboratory professionals follow strict quality procedures to ensure the integrity of your results.



Sample QC

LIBRARY CONSTRUCTION

Library QC

**SEQUENCING** 

Sequencing QC

**RAW DATA OUTPUT** 

Data QC

**BIOINFORMATICS ANALYSIS** 

Delivery QC



# BGI has extensive experience for metagenomics projects

# **Sequencing Service Specification**

DNBSEQ Metagenomics Survey Sequencing services are executed with the DNBSEQ sequencing system.



### Sequencing Service

- 150bp Paired-End sequencing
- · Data output, standard and custom data analysis
- Available data storage and bioinformatics applications



# Sequencing Quality Standard

Guaranteed ≥ 80% of bases with quality score ≥Q30



#### **Turnaround Time**

- Typical 18 working days from sample QC acceptance to filtered raw data availability
- Expedited services are available, contact your local BGI specialist for details








# **DNBSEQ Sequencing System**

DNBSEQ is an innovative high-throughput sequencing solution, developed by BGI's Complete Genomics subsidiary in Silicon Valley. The system is powered by combinatorial Probe-Anchor Synthesis (cPAS), linear isothermal Rolling-Circle Replication and DNA Nanoballs (DNB™) technology, followed by high-resolution digital imaging.

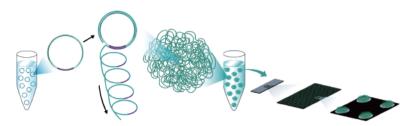

The combination of linear amplification and DNB technology reduces the error rate while enhancing the signal. The size of the DNB is controlled in such a way that only one DNB is bound per active site on the DNBSEQ flow cell. This densely patterned array technology provides optimal sequencing accuracy and increases flow cell utilization.

#### STANDARD ANALYSIS

- · Data filtering
- Alignment
- · Metagenomic De Novo assembly
- Non-redundant gene catalogue
- Prophage transposable element prediction
- Functional annotation based on KEGG, CAZy, eggNOG, CARD
- Species composition and diversity analysis
- Quantitative and differential analysis of gene abundance
- Principal component analysis

#### **CUSTOMIZED ANALYSIS**

Further customization of bioinformatics analysis to suit your unique project is available. Please contact your BGI technical representative.

#### SAMPLE REQUIREMENTS

We can process your extracted genomic DNA from a variety of environmental samples with the following general requirements:

SAMPL	E DNA AMOUNT AND CONCENTRATION	MINIMUM SAMPLE VOLUME
Regula	Intact genomic DNA ≥ 200 ng concentration ≥ 8 ng/µL	15 μL

### **Request for Information or Quotation**

Contact your BGI account representative for the most affordable rates in the industry and to discuss how we can meet your specific project requirements or for expert advice on experiment design, from sample to bioinformatics.

info@bgi.com www.bgi.com

# **BGI Offices**

BGI Americas
One Broadway, 14th Floor
Cambridge, MA 02142,
USA

**BGI Europe** 

Jutrzenki 12 A, 02-230 Warszawa, Poland **BGI** Asia

Building NO.7, BGI Park, Yantian District Shenzhen, Guangdong Province, China

#### **BGI Australia**

L6, CBCRC, 300 Herston Road, Herston, Brisbane, Queensland 4006. Australia

For Research Use Only. Not for use in diagnostic procedures (except as specifically noted).

Copyright© BGI 2022. All trademarks are the property of BGI, or their respective owners. This material contains information on products which is targeted to a wide range of audiences and could contain product details or information otherwise not accessible or valid in your country. Please be aware that we do not take any responsibility for accessing such information which may not comply with any legal process, regulation, registration or usage in the country of your origin. Note, BGI's genetic testing products have not been cleared or approved by the US FDA and are not available in the USA. For Research Use Only. Unless otherwise informed, all sequencers and sequencing reagents are not available in Germany, USA, Spain, UK, Hong Kong, Sweden, Belgium or Italy. Certain sequencing services are not available in USA and Hong Kong. Please contact a representative for regional availability. The company reserves the right of final interpretation.





BGI Genomics

BGI\_Genomics

We Sequence, You Discover