

# 16S/18S/ITS Amplicon Sequencing

## Service Description

Over 99% of natural microorganisms can not be isolated and cultured clonally, while traditional isolation and culture-dependent methods challenge the study of microorganisms in their natural environment.

Metagenomic studies of genetic material that is directly recovered from environmental samples, have benefitted greatly from advanced NGS technology as a method for the exploration of microbial biodiversity.

16S and 18S rDNA are hypervariable regions in the 16S or 18S rRNA genes in bacteria and fungi, while ITS (Internal Transcribed Spacer) is the spacer DNA between the small-subunit and large-subunit rRNA genes in bacteria, fungi and archaea.

Sequence comparison of 16S/18S/ITS regions is widely used in taxonomy and molecular phylogeny because of the easy amplification by PCR, even from low quantities of DNA, while they have a high degree of variation even between closely related species.

## Sequencing Service Specification

BGI 16S/18S/ITS Amplicon Sequencing services are executed with DNBSEQ sequencing system.



### Sample Preparation and Services

- PCR will be used to isolate different 16S/18S/ITS regions
- PE300 sequencing by DNBSEQ
- Sequencing data are available in standard file formats
- Custom bioinformatic data analysis is available
- Data storage services are available



### Sequencing Quality Standard

- Up to  $\geq 80\%$  of bases with a  $\geq 30$  quality score, depending on the chosen sequencing strategy
- Recommended sequencing coverage is dependent on the complexity of the sample




### Turn Around Time

- Typical 35 working days from sample QC acceptance to data availability
- Expedited services are available, contact your local BGI specialist for details

## Project Workflow

We care for your samples from the start through to the result reporting. Highly experienced laboratory professionals follow strict quality procedures to ensure the integrity of your results.

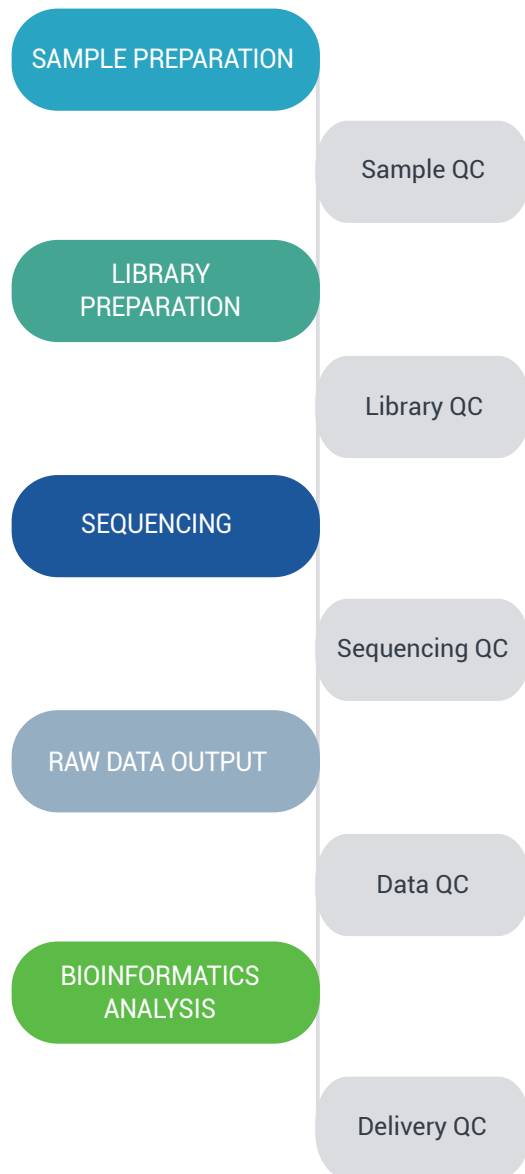

## Data Analysis

Besides clean data output, BGI offers a range of standard and customized bioinformatics pipelines for your sequencing project.

Reports and output data files are delivered in industry standard file formats: FASTQ, .xls and .png

### STANDARD ANALYSIS

- Data filtration
- Overlap paired-end reads to form tags
- Tags are clustered into OTU. PCA, Venn diagram. A rank curve will be generated based on OTU abundance
- Species are classified by OTU annotation, based on which a species profiling histogram, heat map and a phylogeny tree will be provided
- Alpha diversity indexes are generated for in single sample
- Beta diversity and clustering analysis are performed for multiple samples
- Comparative analysis is used to screen significant differences in multiple samples

### CUSTOMIZED ANALYSIS

Further customization of Bioinformatics analysis to suit your unique project is available:

Please contact your BGI technical representative.

## Sample Requirements

We can process intact genomic DNA or PCR fragment samples

	SAMPLE REQUIREMENT	CONCENTRATION
Genomic DNA	≥50ng	≥6ng/μL
PCR Fragment	≥3μg	≥30ng/μL

For Research Use Only. Not for use in diagnostic procedures (except as specifically noted).

Copyright© BGI 2022. All trademarks are the property of BGI, or their respective owners. This material contains information on products which is targeted to a wide range of audiences and could contain product details or information otherwise not accessible or valid in your country. Please be aware that we do not take any responsibility for accessing such information which may not comply with any legal process, regulation, registration or usage in the country of your origin. Note, BGI's genetic testing products have not been cleared or approved by the US FDA and are not available in the USA. For Research Use Only. Unless otherwise informed, all sequencers and sequencing reagents are not available in Germany, USA, Spain, UK, Hong Kong, Sweden, Belgium or Italy. Certain sequencing services are not available in USA and Hong Kong. Please contact a representative for regional availability. The company reserves the right of final interpretation.

## Request Information or Quotation

Contact your BGI account representative for the most affordable rates in the industry, to discuss how we can meet your specific project requirements or for expert advice on experiment design, from sample to bioinformatics.

[info@bgi.com](mailto:info@bgi.com)

## BGI Offices

### BGI Americas

One Broadway, 14th Floor  
Cambridge, MA 02142,  
USA

### BGI Europe

Jutrzenki 12 A,  
02-230 Warszawa,  
Poland

### BGI Asia

Building NO.7, BGI Park,  
Yantian District Shenzhen,  
Guangdong Province, China

### BGI Australia

L6, CBCRC, 300 Herston  
Road, Herston, Brisbane,  
Queensland 4006, Australia



BGI Genomics



BGI\_Genomics



**We Sequence, You Discover**