


Service Description

Protein identification analysis mainly includes gel spot identification and gel band identification, which can be used when it is necessary to identify a specific range of proteins, or when there are only a few kinds of proteins in a sample.

Gel spot identification is mainly used for MS identification of the designated gel spot in two-dimensional (2D) electrophoresis gel. Firstly, the gel spot sample is decolorized and enzymolysed, and then detected by LC-MS/MS. Using the obtained mass spectra, the target protein can then be identified via cross referencing established databases.

Gel band identification is mainly used for MS identification of a protein solution in which protein type is relatively small (dozens to 100) or the designated band in sodium dodecyl sulfate polyacrylamide gel electrophoresis (SDS-PAGE). Firstly, the gel band or protein solution samples are enzymolysed, and then detected by LC-MS/MS. Using the obtained mass spectra, the proteins are then identified via cross referencing established databases.

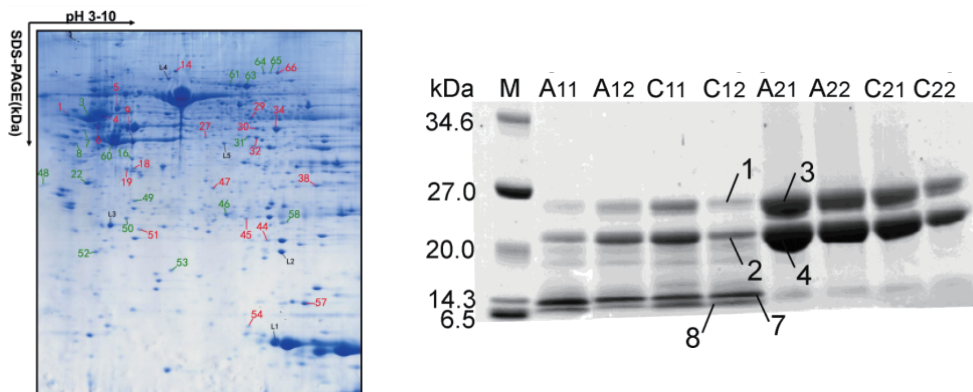

Gel spot (left) and gel band (right)

Zhan X Q, et al., How many proteins can be identified in a 2DE gel spot within an analysis of a complex human cancer tissue proteome, (2018)


BGI has extensive experience in the field of Protein Identification with reliable workflows using nano LC-MS/MS and a bioinformatics infrastructure that is second to none.

Research Applications


- Identification of target proteins
- Identification of protein mixture
- Protein-protein interaction studies

Protein Identification Analysis Workflow

Gel spot identification:

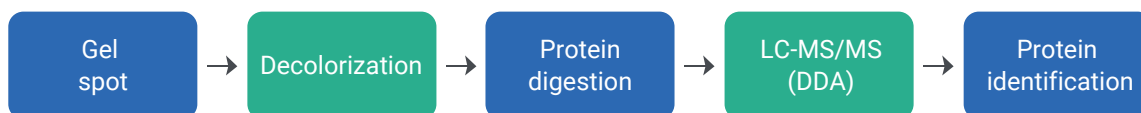


Gel band identification:

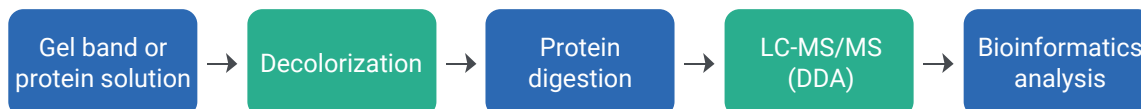

Bioinformatics Analysis Workflow


Standard:

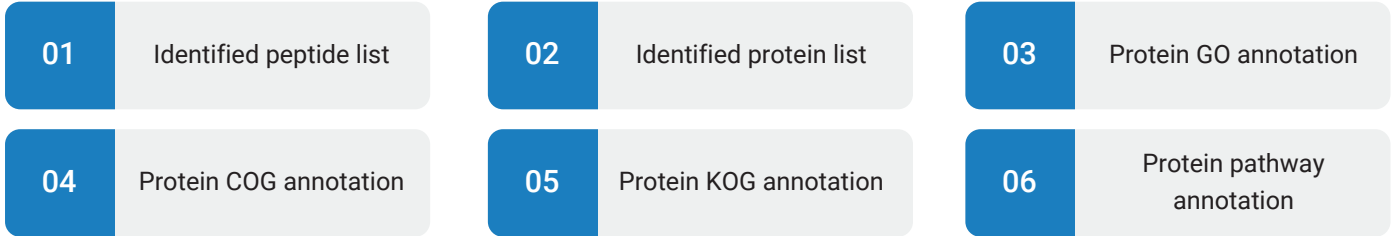


Customized Solutions:

Quantitative analysis of protein relative abundance in a single sample (iBAQ)


Examples of Protein Function Annotation

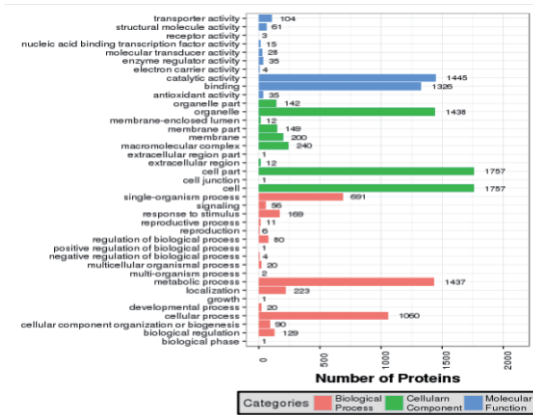


Protein GO Annotation

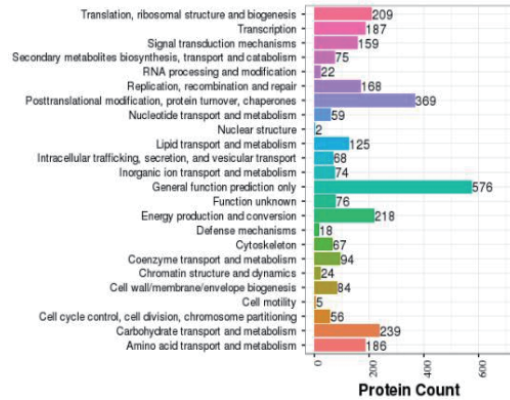

Protein COG Annotation

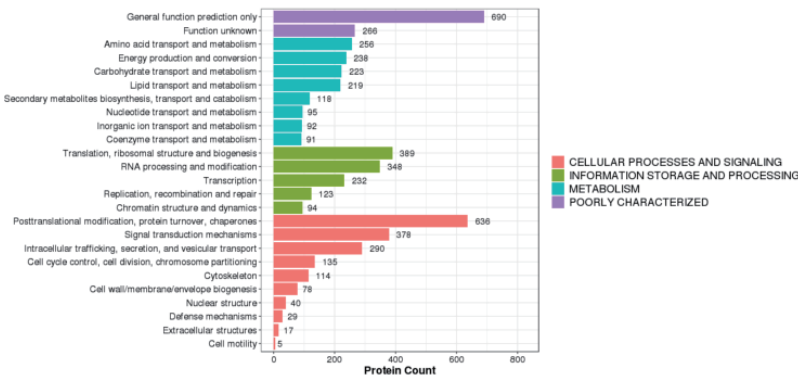

Protein KOG Annotation

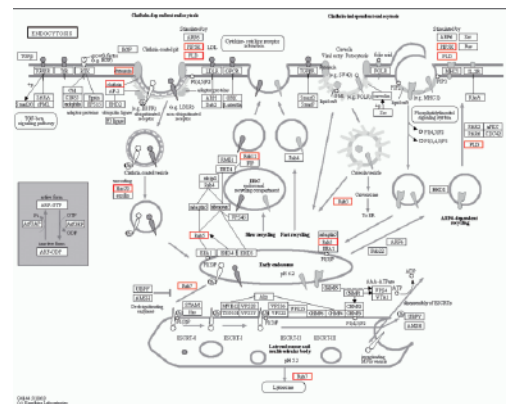

Protein pathway annotation

General Sample Requirements

SAMPLE \ PRODUCT	GEL SPOT IDENTIFICATION	GEL BAND IDENTIFICATION	MODIFICATION SITE IDENTIFICATION
Coomassive or silver stained gel spot/band	Visible spot	Protein $\geq 1 \mu\text{g}$	Protein $\geq 1 \mu\text{g}$. Projects that need enrichment is not accepted.
Single protein (protein solution purified after IP or CO-IP)	/	$\geq 5 \mu\text{g}$, $\geq 0.5 \mu\text{g}/\mu\text{L}$	$\geq 10 \mu\text{g}$, $\geq 0.5 \mu\text{g}/\mu\text{L}$. If enrichment is required, the amount of sample sent needs to be increased by 3 times.

To Learn More

To learn how your research can benefit from BGI's extensive experience in Protein Identification, visit www.bgi.com, write to us via info@bgi.com or contact your local BGI office.

International Head Offices

BGI Americas

One Broadway, 3rd Floor
Cambridge, MA 02142,
USA
Tel: +1 617 500-2741

BGI Europe

Ole Maaløes Vej 3,
DK-2200 Copenhagen N,
Denmark
Tel: +45 7026 0806

BGI Asia

16 Dai Fu Street,
Tai Po Industrial Estate,
New Territories, Hong Kong
Tel: +852 36103510

BGI Australia

L6, CBCRC, 300 Herston
Road, Herston, Brisbane,
Queensland 4006, Australia
Tel: +61 (07) 3362 0475



BGI Genomics



BGI_Genomics

Services For Research Use Only. Not For Use in Diagnostic Procedures (unless specifically noted).

Copyright ©2021 BGI. The BGI logo is a trademark of BGI. All rights reserved.

All brand and product names are trademarks or registered trademarks of their respective holders.

Information, descriptions and specifications in this publication are subject to change without notice.

DNBSEQ™ is a trademark of MGI Co. Ltd.

DNBSEQ™ technology-based services are executed at our service laboratories in China.

Published: February 2021

Version:1.0

